

KemSure[®]

Formaldehyde Neutralizing Powder

Catalog #9144

Directions:

1. Pour KemSure™ Formaldehyde Neutralizing Powder into container with waste Formaldehyde solution which you intend to neutralize (one container neutralizing powder per ¼ gallon waste solution).
2. Stir solution to mix.
3. When reaction is complete (30 minutes), the neutralized solution is then safe for disposal.

Use the entire contents of this container in one application. This product is not designed to be used in smaller portions.

Note: Always maintain adequate ventilation in accordance with applicable regulations. Disposal is subject to federal, state, and municipal sewer regulations.

Precaution:

This product is not to be ingested or brought into contact with the skin and eyes. Review manufacturer's Safety Data Sheet. Keep away from children.