

# *KemSure*<sup>®</sup>

## Formaldehyde Neutralizing Powder

### Catalog #9144

**Directions:**

1. Pour KemSure™ Formaldehyde Neutralizing Powder into container with waste Formaldehyde solution which you intend to neutralize (one container neutralizing powder per ¼ gallon waste solution).
2. Stir solution to mix.
3. When reaction is complete (30 minutes), the neutralized solution is then safe for disposal.

*Use the entire contents of this container in one application. This product is not designed to be used in smaller portions.*

*Note:* Always maintain adequate ventilation in accordance with applicable regulations. Disposal is subject to federal, state, and municipal sewer regulations.

**Precaution:**

This product is not to be ingested or brought into contact with the skin and eyes. Review manufacturer's Safety Data Sheet. Keep away from children.