

Cat. No.
**9520- (P/pail
or T/tote)**
One Kit

KemSure

FORMALDEHYDE NEUTRALIZING SPILL CONTROL KIT

For Use on Formaldehyde Disinfectant Spills up to ½ gallon.

Contents:

- **Two (2) Formaldehyde Neutralizing Spill Response Chemsorb™ Pillows*****
***** Additional Pillows (set of 2) may be purchased with Catalog #9520-S*****
- **One (1) pair of Nitrile Gloves**
- **One (1) AAMI Level 3 Isolation Gown**
- **One (1) Procedure Mask with elastic ear loops**
- **One (1) Full Length Face Shield**
- **One (1) pair Non-Skid Knee-High Boot Covers**
- **One (1) Lightweight Bouffant Cap**
- **One (1) Temporary Disposal Bag with Label**

DIRECTIONS:

1. Open your KemSure® Formaldehyde Neutralizing Spill Control Kit.
2. Remove all contents from kit.
3. Put on gown, booties, gloves, mask, cap, and face shield as necessary.
4. Place one or more Formaldehyde Neutralizing Spill Pillow(s) on spill area. Start at outer edge of spill, apply light pressure and sweep in to absorb all the liquid.
5. Place used pillow(s) in temporary spill bags and seal.
6. Contents should be labeled and management notified. Dispose of material in accordance with Federal, State, and local regulations. **DO NOT REUSE PILLOWS.**
7. Replace Spill Kit -- Re-Order -- Cat. # **9520- (P/pail or T/tote)**. To purchase individual components of this kit, please see the Inventory Card included inside.

DO NOT REUSE PILLOWS

The **KemSure**® Formaldehyde Neutralizing Spill Control Kit is designed for small spills up to ½ gallon. Each Pillow will pick-up approximately 32oz of liquid. Spills should be cleaned up by trained personnel only. Dispose of spilled materials in accordance with Federal, State and local regulations. Avoid generating dust. In case of eye contact, flush with large quantities of water. Read Safety Data Sheet (SDS) for information on product safety and handling procedures.

Chemsorb™ is a Trademark of Precision Laboratories, Inc.

Caution:
**Do not use Chemsorb™ Pillows
for Hydrofluoric acid.**
Keep out of Reach of Children.