

# KemSure™ **FORMALDEHYDE SPILL CONTROL KIT**

**Provides safe and easy clean up of formaldehyde spills**

## CONTAINS

- 2 - Formaldehyde Neutralizing Spill Response Chemsorb Pillows
- 2 - Temporary Spill Disposal Bags
- 1 - Pair Nitrile Chemical Resistant Gloves
- 2 - Identification Labels.

**Lot #** F010

**Expiration Date:** Dec 2012



Telephone: (631) 454-6565 • (800) 553-0330  
Fax (631) 454-8083

## CAUTION

**Do not use Chemsorb for hydrofluoric acid.**  
*See Directions for other precautions*

**Catalog Number 9145**

One Kit

Made in USA

## **FORMALDEHYDE SPILL CONTROL KIT**

Each FORMALDEHYDE SPILL Pillow will neutralize approximately ¼ gallon of formaldehyde liquid.

**Spills should be cleaned up by trained personnel only.**

**Dispose of spilled materials in accordance with federal, state and local regulations.**

### *Directions For Use*

1. Open your KemSure™ Formaldehyde Spill Control Kit. It should contain:

2 - Formaldehyde Neutralizing Spill Response Pillows	2 - Temporary Spill Disposal Bags
1 - Pair Nitrile Chemical Resistant Gloves	2 - Identification Labels.
2. Remove all contents from kit. Put on gloves.
3. Place one or more Formaldehyde Neutralizing Spill Pillow(s) on spill area. Start at outer edge of spill, apply light pressure and sweep in to absorb all the liquid.
4. Place used pillow(s) in temporary spill bags and seal.
5. Contents should be labeled and management notified. Dispose of material in an appropriate and responsible manner. **DO NOT REUSE PILLOWS.**
6. **Replace** Spill Kit, -- re-order -- *Cat. # 9145*

*Kem Medical Products Corp.*

Telephone: (631) 454-6565 • (800) 553-0330  
Fax: (631) 454-8083

## CAUTION: KEEP OUT OF REACH OF CHILDREN.

**Do not use Chemsorb for hydrofluoric acid. Avoid generating dust. In case of eye contact, flush with large quantities of water. Read Material Safety Data Sheet (MSDS) for information on product safety and handling procedures.**

CHEMSORB® is a Registered Trademark of Precision Laboratories, Inc.